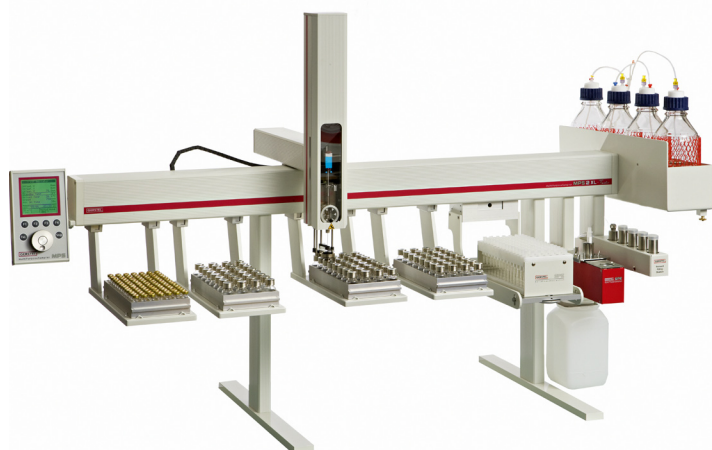



**GERSTEL**


## Solid Phase Extraction

# SPE

## Specifications

### Uses

Sample preparation and sample introduction system for automated Solid Phase Extraction SPE based on standard cartridges. Conditioning, extraction and elution is performed in a sealed system. An automated concentration can be selected for improved detection limits. SPE can be combined with introduction of the eluate to a GC or LC system or performed off-line for subsequent analysis of the SPE eluate using a separate GC/MS or LC/MS system.

### System configuration

- compatible with most standard GC and LC systems
- automated operation based on the GERSTEL MultiPurpose Sampler MPS
- GERSTEL Solvent Filling Station SFS included

### Automated SPE steps

- cartridge conditioning
- sample loading
- cartridge drying
- elution
- eluate evaporation and analyte concentration
- sample reconstitution
- solvent exchange

### Cartridge types

- standard SPE cartridges with GERSTEL transport adapter
- 1 mL, 3 mL and 6 mL
- height of packing inclusive of frit  
4.5 ... 23 mm (for 1 mL and 3 mL cartridges)  
9.5 ... 23 mm (for 6 mL cartridges)
- more detailed information is available in a separate flyer

### Capacity

- 72 × 6 mL cartridges
- 98 × 3 mL cartridges
- 98 × 1 mL cartridges

### Sample vials

- 2 mL, 4 mL or 10 mL

### Sample volume

- max. 8 mL per vial
- a sample can be split up into several vials

### Eluate collection vials

- 2 mL, 4 mL or 10 mL

### Syringes

- 1 mL or 2.5 mL, with gas connection
- syringe volume does not limit the injection volume, multiple injections are possible

### Injection volume

- 10% or more of total syringe volume
- in combination with LC, smaller sample volumes can be introduced using injection valves with small sample loops

### Loading speed


- 10 ... 250 µL/s

### Evaporation temperature

- ambient ... 120 °C

### Agitation speed

- 250 ... 750 rpm

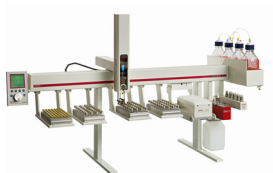

## Solid Phase Extraction SPE

### Control

- in combination with the GERSTEL MAESTRO software, integrated in the Agilent® Technologies ChemStation software or operated in stand-alone mode
- only one method and one sequence table required for the complete system including GC/MS or LC/MS system when integrated in the ChemStation software
- sample preparation steps can be selected by mouseclick from a drop-down menu
- SPE and chromatographic analysis of the preceding sample are performed simultaneously ensuring optimal productivity and throughput

### Operating conditions

- 15 ... 35 °C
- relative humidity max. 50-60%, non-condensing
- max. 4615 m above sea level

### Storage conditions

- -20 ... 50 °C
- relative humidity max. 50-60%, non-condensing
- max. 4615 m above sea level

### SPE Station

- dimensions 22.5 x 9 x 26 cm (H x W x D)
- weight 1.7 kg