




Soprane II


License / Trial activation



The first thing to do is to download the software, you can download it with the following link: <https://www.srainstruments.com/p/soprane-ii/>, click on “Trial version” (it’s the complete installation setup).


Then execute the setup installation file.

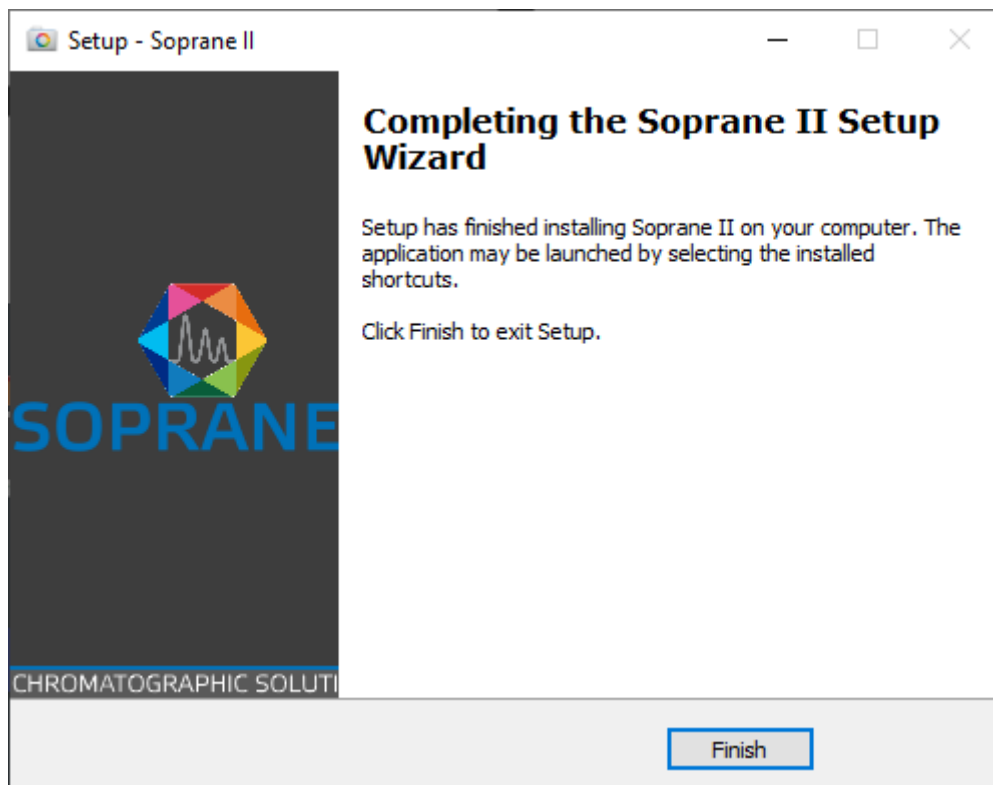

Once *Soprane II* is correctly installed, you will see a new icon on your desktop  this is the configurator, you can also find it in the installation folder under the name *SRA.Soprane.Configuration.exe*.


Start the configuration application, wait a few seconds and the message “*Invalid license*” confirms that the license is not validated.



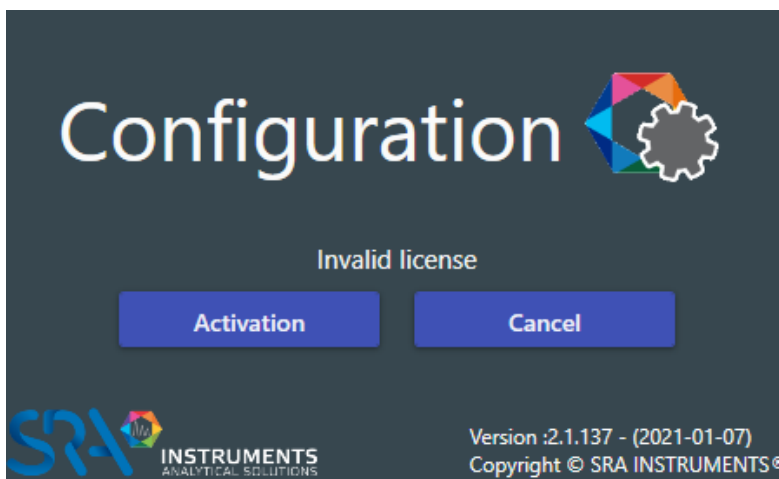


Click on the **Activation** button; the following window appears.

If the configuration starts without showing the activation button, please close the configuration and open the following application: "C:\Soprane \SRA.Soprane.LicenseActivation.exe"

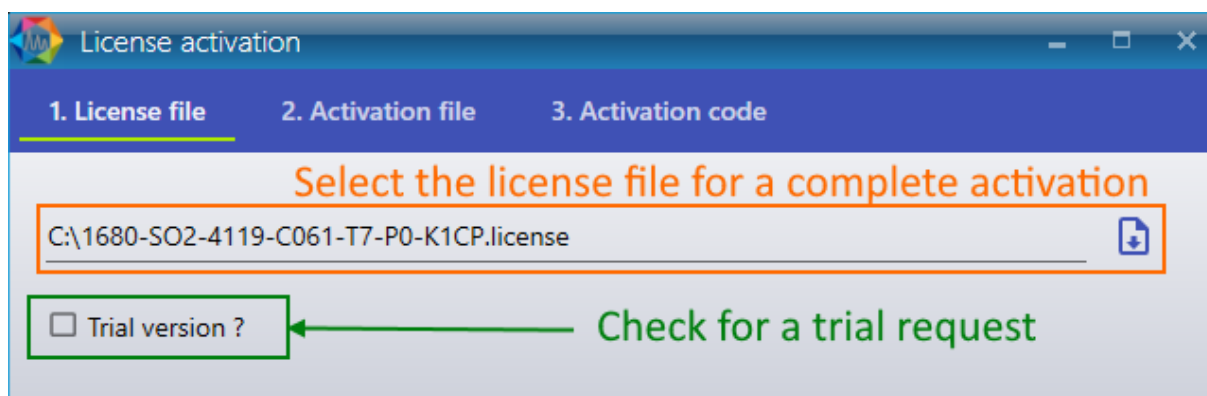


- For a license activation, select the "license" file sent by SRA Instrument.
- For a trial activation just check "Trial version", and leave the license field empty.

Once one of these field is filled, the next tab will be automatically selected.

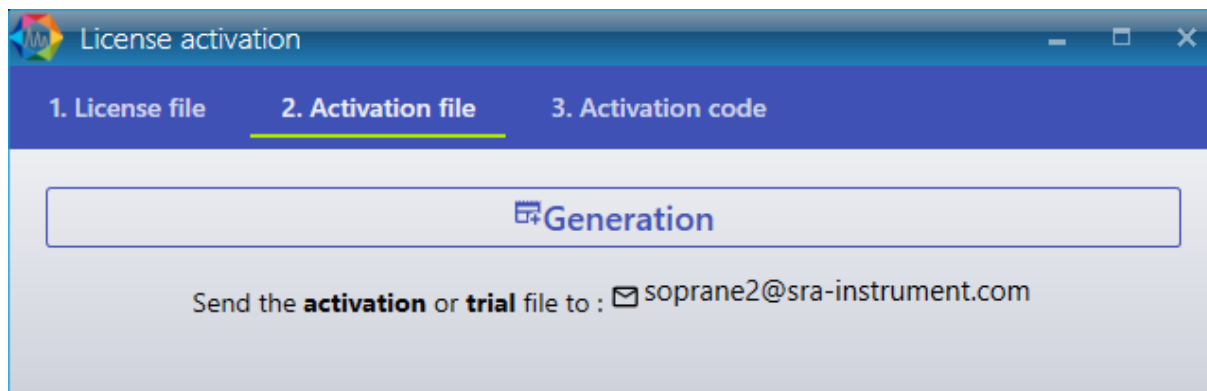

Click on **Generation**. A file will be created in Soprane’s folder with the extension “.skey”. Send this file by e-mail to soprane2@sra-instruments.com.

SRA Instruments will send you the activation code you need. Then fill in the field **Validation** with this code (example a6d457c2-2f61-89e2-6833-eed960c74897).

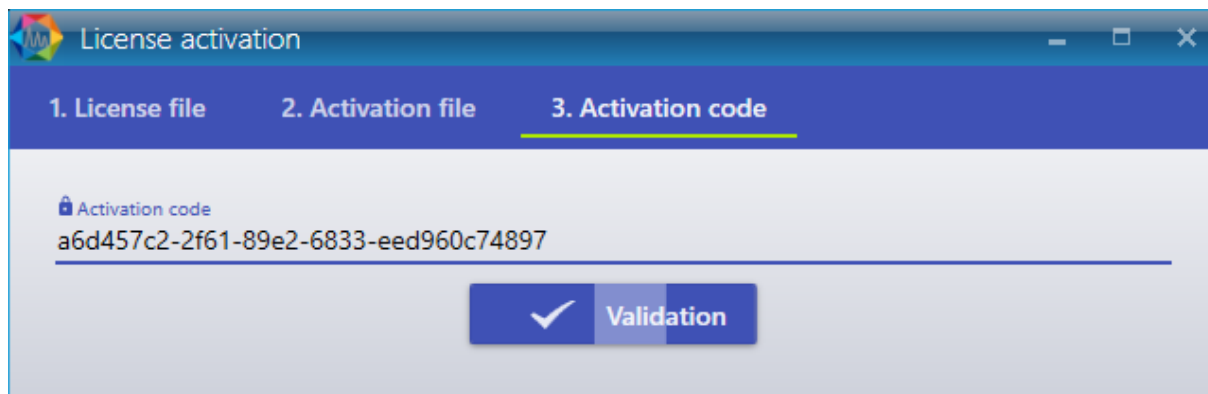

Click on Validation, if the code is correct, your 30 days trial or your license will be activated.

Now enjoy all the new features of Soprane II!

